


## Dilucup Elegance NaCl

**Product No.** 40DENACL3 (2x 16 x 21 cups), 40DENACL6 (2 x 8 x 42 cups)

**Use and description:** Ready-to-use plastic cups, each contains 9 mL medium to be used for diluting microbiological samples together with a special shaker, Dilushaker. By adding a sample of 1 mL to the first cup a dilution of 1:10 will be achieved. 1 mL from this cup is then transferred to the next to get a dilution of 1:100. The procedure is continued until the required dilution is reached

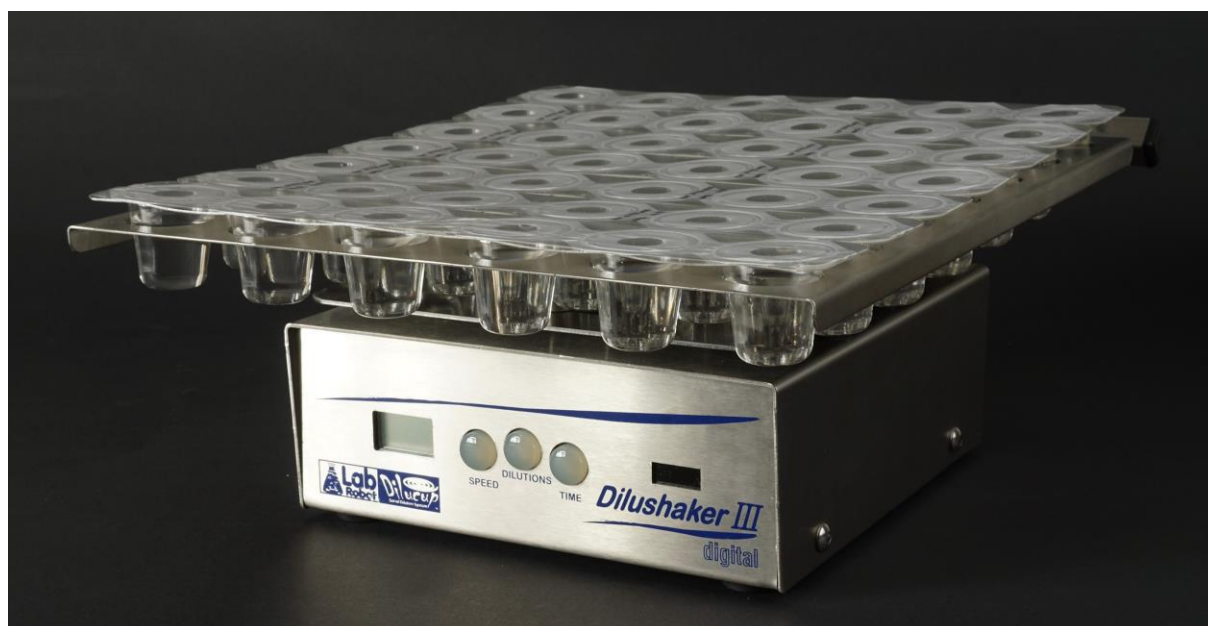

**Composition per litre:**

NaCl	8,5 g
Buffers	1,4 g

**pH:** 7,0 +/-0,2 at 25 °C

**Volume per cup:** 9 mL +/- 2 %

**Sterilization:** E-Beam > 25kGy

**Material of the cups:** PET (recyclable)

**Expiry date:** When unopened >18 months after production

**Storage:** Unopened: +2 to +25°C. +2 to +8°C when the box is open

**Packaging:** The cups are delivered in blisters of 7 pre-cut rows of 3 cups (40DENACL3) or 6 cups (40DENACL6). Each box contains 672 cups distributed in 2 bags of 336 cups

**Weight and volume:** 8,5 kg/box - 0,032 m<sup>3</sup>/box